



- 1) ASSEMBLE THE 4" FRONT AND BACK RAILS, USING STUD PLATES AND LOCK NUTS PROVIDED. BEFORE TIGHTENING NUTS, MAKE SURE THE FRONT AND BACK RAIL ARE FLUSH TO THE TOP OF THE LEG ASSEMBLIES AND PUSHED INWARD APPROXIMATELY 1/8" TO ALLOW THE WORKSURFACE TO BE MOUNTED PROPERLY. NOTE: HOLES IN RAILS SHOULD BE FACING UP.
- 2) CLEAN THE ENTIRE UNDERSIDE OF THE WORKSURFACE AND THE TOP RAILS OF THE FRAME WELL USING RUBBING ALCOHOL.
- 3) PUT A BEAD OF LIQUID NAILS ALONG THE TOP RAILS OF THE ENTIRE FRAME.
- 4) CAREFULLY PLACE THE EPOXY RESIN TOP FLUSH WITH THE BACK LEGS AND EVEN ON THE SIDES. LET LIQUID NAILS DRY OVERNIGHT

PROLINE  
EPOXY RESIN HDL BENCH ASSEMBLY  
DRAWING # A7650 - ER